

1. Blackground, Darkfield microscope, the biggest expectation of customers




- ① Base Assembly include: **CCD color camera and microscope.**
- ② 2 Ocular , 4 Object lens: 4X、 10X、 40X、 100X 4x and 10x is for the bright background, 40x and 100 x for the Black ground.

② In general, darkfield microscope need an additional darkfield condenser to complete its function. When the operator tests live blood, has to set the darkfield condenser first. After that, he has to take off the darkfield one, uses the brightfield condenser to observe dry blood. The operational procedure is complex.


Our microscope, has obtained three patents, not only display a very clearly image under the darkfield, but also freely switch between darkfield and brightfield.

There is only one adjustor ring under the conserder of our Blackfield microscope. The adjustor ring puts in the Position A when starts to observe the live blood. We simply turn the adjustor ring to Position B then can test dry blood under the brightfield. The whole procedure is very easy. There is no need to purchase other device and the improvement decreases your cost and saves your time.

Clear image, simple operation, low cost, aren't those the biggest expectations of our customers?

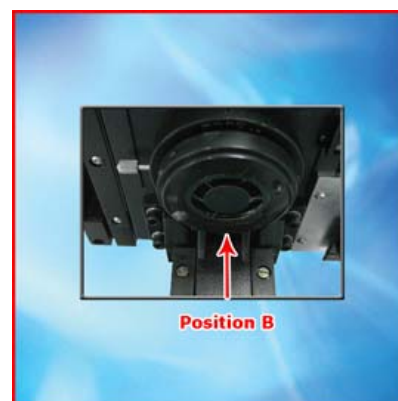

2. Galleries

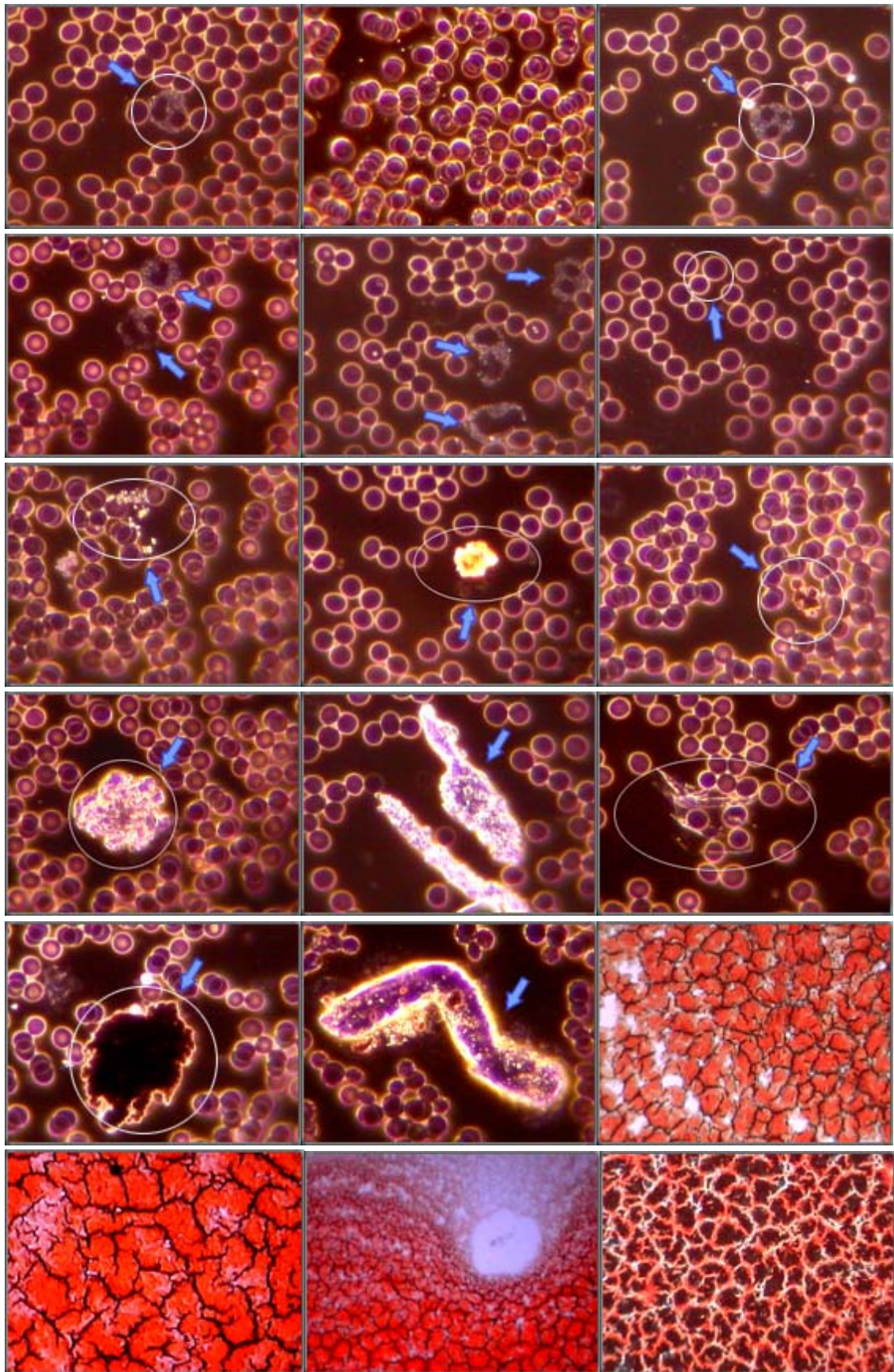

3. Free Accessories include:

- ① **Aluminum alloy carrying case with foam liner; (570x270x400mm 20kg)**




- ② **External capture card;**



- ③ **Software and software dog;**



- ④ **Box of 50 blank microscope slides;**

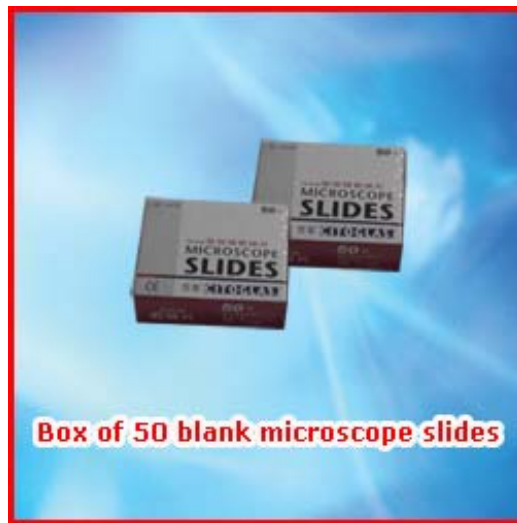

- ⑤ **Bag of 100 lancets;**



- ⑥ **Lancing device;**



⑦ 100 pieces cover slips;



⑧ Rubber bulb;



⑨ Halogen bulb (6V,20 Watts) and Tube fuse.;



⑩ Lens paper;



⑪ Training book

Catalogue	
● Illustration	1
● Terminology	2
● Operation	6
● Additional Points	13
● Contact us	17

Read this manual before using the microscope

⑫ Dust cover;